testoval[®]

NITRIT-TEST

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Directions for Use

Remove the test cell from the comparator and fill it up to the mark with water sample. Add 2 dosing spoons full of nitrite reagent and dissolve by stirring with the red spoon. After 15 minutes, reinsert the test cell into the comparator, hold the comparator against a bright background or light source, and match the colour intensity of the central field (test solution) with one of the colours in the peripheral colour fields. Read the value in ppm NO₂, which is printed under the matching colour field.

Should the colour in the central field be more intensive than any of the colours in the peripheral fields, repeat the test with a water sample diluted 1:10. For this purpose fill the graduated plastic tube up to the lower mark with sample water and dilute with nitrite-free water (e. g. distilled water) to the upper mark. Transfer the diluted sample into the test cell, which has been rinsed with nitrite-free water and proceed as described above. The value obtained must in this case be multiplied by 10.

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